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RUBBER STOPPER/ ALUMINUM CAP WASHING MACHINE  
Professional pharmaceutical equipment and overall solution supplier

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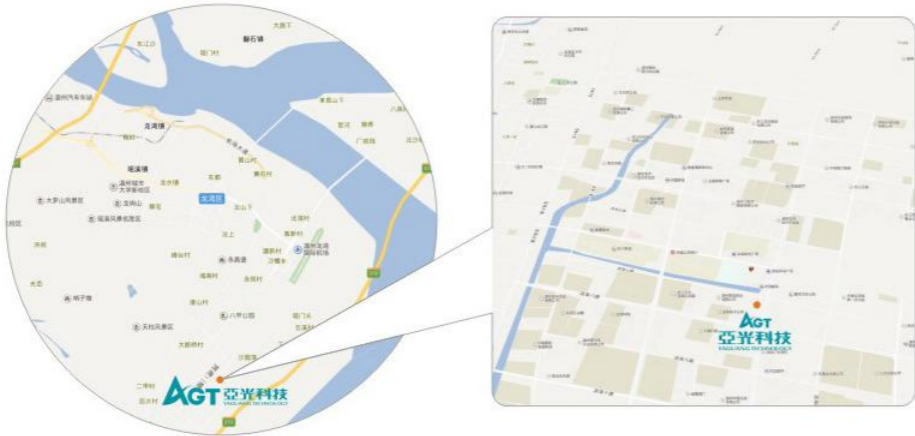
English website

亚光工业园 Yaguang Industrial Zone

Zhejiang Yaguang Technology Co., Ltd is a national certified high-tech enterprise specializing in the design, production and sales of pharmaceutical equipment since established on 1996. Yaguang has been focusing on pharmaceutical equipment industry for years and developed the first Rubber Stopper Washing Machine in 1997, the first Agitated Nutsche Filter Dryer and other core equipment of pharmaceutical process system in 2000, Yaguang has a leading position in China's pharmaceutical industry.

The main products of Yaguang fall into three categories: Washing Machine (including Rubber Stopper / Aluminum Cap Washing Machine), API production line (Reaction Tank, Crystallizing Tank, Agitated Nutsche Filter Dryer, Conical Pulse Mixing Dryer, Split Charging Weighing System and so on), Efficient Rotating Distillation Bed. More than that, Yaguang is the drafting unit for industry standards of Rubber Stopper Washing Machine and Agitated Nutsche Filter Dryer, has obtained more than 100 patented technologies, the equipment has passed certifications such as American FDA, EU cGMP, Japanese GMP in a number of domestic and abroad pharmaceutical industries.

Yaguang has a provincial R&D center with manufacturing licenses of D1, D2 pressure vessel in China. With the solid technical foundation, advanced management concepts and excellent corporate culture, Yaguang is developing with a steady pace and being a global professional pharmaceutical process program supplier.

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Advanced technology and excellent quality facilitate the development of pharmaceutical and chemical equipment in China

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Our company always regards talent cultivation as the foundation for enterprise development. Through trainings at various levels, in multiple forms and from varied channels, we cultivate all kinds of talents. Yaguang continuously introduces experienced senior and intermediate talents, enhances talent storing, reinforces combination of theory and practical experience and provides customers with services of "sincerity, fairness, high quality and high efficiency" with our powerful technical strength.

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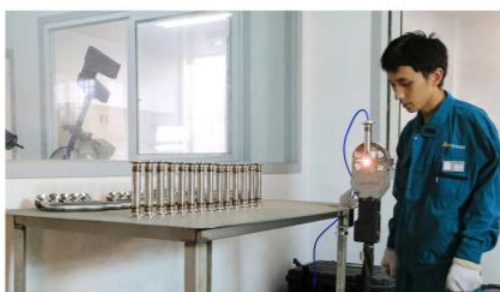
The basis of market competition, inseparable from the production capability, and can not do without an advanced hardware facilities. Fully aware of the importance of hardware facilities, Yaguang takes the lead in the introduction of foreign advanced production equipment in the industry; during recent years Yaguang has further invested in updating the hardware facilities, to produce high-end quality products with the internationalized production facilities to meet the GMP standard.

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### Products

1. Washing machine series: The main products include medical rubber stopper washing machines, aluminum cap washing machines and etc. Yaguang is the drafting unit of national industrial standard of rubber stopper /aluminum cap washing machines.
2. Solid-liquid separation and powder treatment series: The main products include Aseptic Dissolving Tanks, Crystallizing Tanks, Agitated Nutsche Filter Dryer, Comminutor, Mixers, Automatic Weighing and Split-charging Systems. Yaguang is the drafting unit of national industrial standard of pharmaceutical Agitated Nutsche Filter Dryer.
3. Solvent recovery: Distillation Bed.

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### Services

Design, manufacturing, installation, commissioning and technical support for the equipment.

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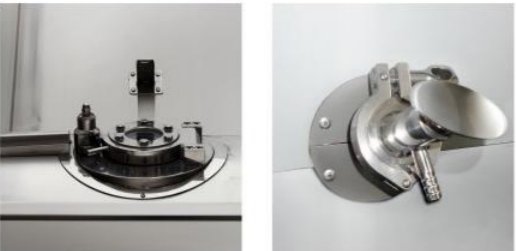
KJ Series Rubber Stopper/Aluminum Cap Washing Machine

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Developed since 1996, the rubber stopper/aluminum cap cleaning, sterilizing and drying equipment has been developed to the sixth generation, forming two major series of 80 specifications. We are the drafting unit of national standards of pharmaceutical rubber stopper/aluminum cap washing machine.

观察/进料口 Observe / feeding nozzle    取样口 Sampling nozzle

— 胶塞清洗机  
Rubber Stopper Washing Machine

进料斗 Feeding hopper

出料口 Discharging nozzle

— 铝盖清洗机  
Aluminum Cap Washing Machine

清洗机种类、功能 Type and Function

种类 Type	适用产品 Applicable Products	主要功能 Main Functions
KJCS超声波胶塞清洗机 KJCS Ultrasonic Rubber Stopper Washing Machine	粉针胶塞、冻干胶塞、水针西林瓶胶塞、大输液胶塞  Powder Injection Rubber Stoppers, Lyophilization Rubber Stoppers, Water Injection Antibiotic Bottle Rubber Stoppers, Large Infusion Rubber Stoppers	自动进料、清洗、取样、硅化、灭菌、干燥、冷却、自动出料、CIP、SIP，全过程电脑自动控制操作  Automatic Feeding, Cleaning, Sampling, Siliconizing, Sterilizing, Drying, Cooling, Automatic Discharging, CIP, SIP, and Whole Process Automatically Computer Controlled.
KJQS气冲式胶塞清洗机 KJQS Air Impact Rubber Stopper Washing Machine		
KJSL铝盖清洗机  KJSL Aluminum Cap Washing Machine	药用铝盖、铝塑复合盖、针管胶塞和注射活塞。  Medical Aluminum Caps, Aluminum-plastic Compound Caps, Needle Tube Rubber Stoppers and Injection Pistons	人工进料、清洗、取样、灭菌、干燥、冷却、自动出料、CIP、SIP，全过程电脑自动控制操作  Manual Feeding, Cleaning, Sampling, Sterilizing, Drying, Cooling, Automatic Discharging, CIP, SIP, and Whole Process Automatically Computer Controlled.

KJ型清洗机基本技术参数 Technical Parameters of KJ series

型号 Model	产量(只) Output (piece)	水箱容积(L) Water Tank Volume	长(mm) Length	宽(mm) Width	高(mm) Height	重量(kg) Weight
KJ-1	10000	330	2200	1650	2200	1800
KJ-2	20000	620	2450	1850	2200	2200
KJ-4	40000	760	2650	1850	2200	2500
KJ-5	50000	925	2850	1850	2200	2500
KJ-6	60000	1130	3300	1850	2200	2700
KJ-8	80000	1380	2850	2100	2300	3000
KJ-10	100000	1630	3050	2100	2300	3200
KJ-12	120000	1930	3350	2100	2300	3400
KJ-15	150000	2280	3100	2450	2500	4000
KJ-18	180000	2630	3350	2450	2500	4800
KJ-20	200000	2940	3700	2450	2500	5000
KJ-22	220000	3350	3650	2600	2600	5200
KJ-24	240000	3450	3750	2600	2600	5300
KJ-26	260000	3650	3950	2600	2600	5500

注：产量中胶塞以Φ19×8.5粉针西林瓶胶塞为标准，Φ20复合铝盖与Φ19粉针胶塞充装率比约为1.5~1.7：1  
Note: Production capacity of rubber stoppers calculates by the powder injection antibiotic bottle of Φ19×8.5, the filling rate between Φ20 composite aluminum cap and Φ19 powder injection rubber stopper is about 1.5~1.7:1.

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### KJ Series Rubber Stopper/Aluminum Cap Washing Machine

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##### Automatic closed discharging device

- ◆ Commonly known as "elephant nose discharging device", can effectively avoid outer environmental pollution when discharging.
- ◆ CIP & SIP.
- ◆ Rubber stopper can be transferred in the confined environment with various docking systems.

##### Central spray and surround 360° spraying technology

- ◆ Central spray is by using high pressure water to spray rubber stopper closely, which is the most effective way to wash pyrogen.
- ◆ Surround 360° spraying technology is applied to equipment self-cleaning. Spray tube rotates along with inner vessel, the spray nozzles spreading on the tube spray and clean barrel and inner vessel to achieve the real no dead corner CIP technology.

##### Patented Back-Door pullout structure

- ◆ Rotating barrel and back-cover totally pull out technology (pioneer in China, patent technology), more meet FDA and EU standard.
- ◆ Thoroughly solve the problem of jacket repair and washing which troubles drug factory for many years.
- ◆ The maintenance work of movement for the rear cap can be processed in the cleaning area of D level (ISO08 level), avoiding from the pollution of the C level cleaning area (ISO7 level) or the B level cleaning area (ISO5 level).
- ◆ The construction will be more reasonable and simplification. It is conforming to the cleaning standard and creating the new ideas of cGMP standard.

##### Electrical control fully meet processing parameters historical record requirements

- ◆ Control system consists of industrial touch panel PC (Panel PC) + Programmable Logic Controller (PLC). It's easy to control full automatically. Operating parameters can adjust in dialog box and process can display dynamically. The alarm system can provide alarm information and carry out parts interlock to prevent wrong operation.
- ◆ Control software use FDA-approved Italian Movicon system. It can carry out rights management, automatic detection of chamber leakage and automatic calibration of temperature and pressure. It can joint to corporate network and easy to improve management and monitoring for equipment and production.

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DQ Series washing machine is non-pressure vessel equipment for the washing and drying of rubber stopper and aluminum cap. The products including DQXT series (with vacuum, electric heating drying function) rubber stopper washing machine, DQXL series aluminum cap washing machine.

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##### Main Features：

Manual feeding, washing, sampling, (siliconizing), drying, cooling, automatic discharging for spiral reversal and CIP.

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
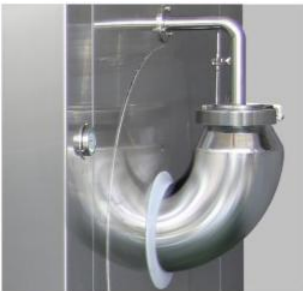
Note: Production capacity of rubber stoppers calculates by the powder injection antibiotic bottle of  $\phi 19 \times 8.5$ ; Production capacity of aluminum caps calculates by the size of  $\phi 20 \times 8.2$

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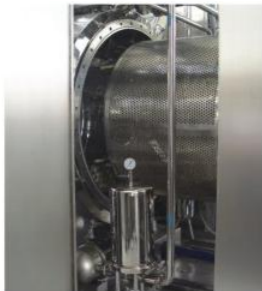Under the protection of isolated orabs laminar flow, there are four process transfer modes according to customer demand

◆Orabs scheme of single aseptic bag automatic packaging and discharging transfer: automatic sealing machine

The main functions include: automatic vacuum bag grabbing, automatic transfer bag placing, automatic bag clamping, automatic vacuum bag opening, discharge metering control, automatic bag sealing, automatic bag discharging, full automatic operation.

◆Orabs scheme of double aseptic bags and single bag packaging and discharging: semi automatic sealing machine

The main functions include: Taking the inner and outer bags in turn manually, operating and starting the outer and inner bags, manually moving the sealing machine back and forth, and automatic heat sealing. The whole process adopts manual and semi-automatic operation.

◆Orabs scheme of double aseptic bag single automatic double package discharge transfer: automatic sealing machine + semi-automatic sealing machine, The main functions include: automatic vacuum grasping inner bag, automatic transfer and placing inner bag, automatic clamping inner bag, automatic vacuum opening inner bag, discharge metering control, automatic inner bag sealing, automatic out inner bag, automatic lifting and transferring sealed inner bag; The outer bag is taken manually, the outer bag is clamped by operation, the inner bag sealed on the lifting plate is put into the outer bag manually, the outer bag is automatically heat sealed, and the outer bag is automatically discharged. The inner bag in contact with materials is operated automatically in the whole process, The manual operation of the outer bag is semi-automatic.

◆Double aseptic bag, double automatic, double encapsulation, discharging and transferring orabs scheme: automatic sealing machine + automatic sealing machine

The main functions include: automatic vacuum grasping inner bag, automatic transfer and placing inner bag, automatic clamping inner bag, automatic vacuum opening inner bag, discharge metering control, automatic inner bag sealing, automatic out inner bag, automatic lifting and transferring sealed inner bag; Automatic vacuum grab outer bag, automatic transfer put outer bag, automatic clip outer bag, automatic vacuum open outer bag, manual put the inner bag sealed on the lifting plate into the outer bag, automatic outer bag sealing, automatic push out the sealed sterile outer bag. Automatic operation is adopted in the whole process of inner and outer bag packaging.

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- ◆The inside and outside aseptic bags are automatically grasped and opened with a fully automatic structure to avoid the risk of manual gloves clamping and opening the bags and polluting the inside bags. In addition, automation can reduce labor intensity.
- ◆The automatic heat sealing machine can move forward and backward – when heat sealing, it is not facing the outlet of the cleaning machine, so as to avoid polluting the rubber plug in the cleaning machine.
- ◆Stepless frequency conversion control of fan in clean unit: always keep the required wind speed in laminar flow protection.
- ◆Glove door misoperation alarm – because the glove door opening in the production process will affect the micro positive pressure in the isolation cabinet and damage the internal class a environment, the glove door is set with misoperation alarm.

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Sterile isolated transfer system for rubber stopper or aluminum cap

## 2 移动层流车转运 Movable laminar flow car transfer

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### System components:

- ◆ Isolation cabinet
- ◆ Transfer barrel
- ◆ Pinch-prevent butterfly valve
- ◆ Movable laminar flow car

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### Solution summary：

Put the barrel in the isolation cabinet by the discharging side. Open the pinch-prevent butterfly valve. Start the discharging to make the rubber stoppers fill in to barrel. After the barrel is full, close the butterfly valve, barrel transferred to laminar flow car. Laminar flow car transfer the barrel to filling machine and link to it .Open the side door of the filling line, lift the material barrel to the same height with feed platform of filling line, transfer the barrel into the oRABs and then move to next process.

### System feature:

- ◆The whole transfer process always under the sterile clean A grade protection condition, avoid exposure pollution
- ◆Isolation cabinet link to laminar flow car without any crack, avoid ambient pollution.
- ◆Roller of the laminar flow car can be lifted, reduced personnel operation when linking to the filling line.

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#### Solution summary：

Use active oRABs, transport rail and transit hopper to transfer rubber stoppers.  
Under the protection of oRABs, motor drive the belt, belt guide the car to finish the hopper transmission, sensitive to the hopper, motor stops rotating, start discharge procedure. After discharge, motor reverse rotate to guide the hopper to the buffer hopper of filling line. Open the barrel door automatically (manually), automatic discharge. Next circulation begins when the discharging is over.

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#### System components:

- ◆ Isolation cabinet
- ◆ Transfer hopper  
(automatic or manual discharging)
- ◆ Transfer car
- ◆ Drive device

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#### System features：

System requirements: Cleaner, filling machine and capping machine produce one to one. Discharging nozzle, filling and capping machine are all in the same line.  
◆ Discharging from the cleaner to the filling line is always under A grade protection. Avoid exposure pollution.  
◆ Whole transmission process is automatic, reduce personnel interference.  
◆ Photoelectric switches control  
◆ Lift feeding speed is adjustable

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#### Solution summary：

It is adopt for toe to toe of rubber stopper washing machine and filling machine. Connect by sealing stainless steel. It used the altitude difference of them. Rubber stoppers drop to filling machine.

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#### Features：

Discharge by permanent discharging structure of elephant nose. The discharging mouth can be wash, sterile, dry and cool individually or with the washing tank together.  
◆ It used the 360° rotation sprayer to wash. There are no dead space since the spray can cover the whole inner face of the transfer pipe.  
◆ The material of the pipe is 316L stainless steel. There are no corner angle and dead space. The polish rate is under  $0.4 \mu m$ .  
◆ It adopts PLC control system. The working process and the time, temperature and pressure can be show on screen.  
◆ It is effective, risk-free and capital intensive, and meets the requirement of new GMP.

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© 亚光企业(AGT INDUSTRY)版权所有